

Compendium of Analytical Laboratory Methods for Food and Feed Safety: Chemical Analytical Manual (CAM)

METHOD NUMBER: C-005.03

POSTING DATE: original posting 2/1/2019; Chemistry Research Coordination Group

posting extension, 12/2020.

POSTING EXPIRATION DATE: 1/31/2023

PROGRAM AREA: Chemical Contaminants and Toxins

METHOD TITLE: Determination of pentobarbital in ingredients of animal origin and in finished pet foods using liquid chromatography tandem mass spectrometry.

VALIDATION STATUS: Level 2, Single laboratory validation

AUTHOR(S): Tara J. Nickel and Christine R. Casey, US FDA Office of Regulatory Affairs/Office of Regulatory Science/Office of Food and Feed Laboratory Operations/ Denver Laboratory

METHOD SUMMARY/SCOPE:

Identification and quantitative determination of pentobarbital in tallow and other grease products of animal origin using Liquid Chromatography-Tandem Mass Spectrometry (LC-MS/MS), using the negative electrospray ionization (ESI) mode. Pentobarbital is extracted from a homogenous portion of tallow, wet dog food, dry dog food, or horsemeat using acetonitrile. After centrifugation, the supernatant is diluted 1:1 with water and analyzed via LC-MS/MS using a solvent standard curve. Deuterated pentobarbital (pentobarbital-D5) is used as an internal standard (I.S.) to correct for sample matrix suppression and/or loss of analyte. Identification of pentobarbital in a sample is based on both correlation of pentobarbital chromatographic retention time (RT) with that of a standard and ion ratio match.

Analytes(s): Pentobarbitol

<u>Matrices:</u> Tallow and other grease products of animal origin, horsemeat, wet (canned) dog food, and dry dog food, wet (canned) and dry cat food, meat/bone meal, and meats including horse, beef, lamb, pork, goat, venison, bison, elk, rabbit, liver, and kidney.

REVISION HISTORY: Revised 7/18/2019. Revision expanded the scope of method from tallow and grease products to include horsemeat, wet (canned) dog food, and dry dog food.

Revised 2/1/2020: Scope was expanded to include wet (canned) and dry cat food; meat/bone meal, and meats including horse, beef, lamb, pork, goat, venison, bison, elk, rabbit, liver, and kidney.

FOOD AND DRUG ADMINISTRATION OFFICE OF REGULATORY AFFAIRS

Denver Laboratory

Document Number: WI-000140

Compendium method C-005.xx

Revision #: 04 Revised: Nov. 2019

Title:

Determination of pentobarbital in ingredients of animal origin and in finished pet foods using liquid chromatography tandem mass spectrometry.

Page 1 of 34

Sections in This Document

1.	Purpo	ose	2
2.	Scop	e	2
3.	Refer	rences	2
4.	Proce	edure	3
	4.1.	Summary of Method	3
	4.2.	Preparation of Samples	4
	4.3.	Preparation of Standards	5
	4.4.	Reagents	8
	4.5.	Equipment (equivalent equipment may be substituted)	8
	4.6.	Instrumentation (equivalent instrumentation may be substituted)	9
Ana	alyte		9
	4.7.	Extraction	11
	4.8.	Data Analysis and Quality Acceptance	12
5.	Gloss	sary	16
6.	Reco	rds	16
7.	Docu	ment History	17
8.	Chan	ge History	17
9.	Attac	hments	17
	Attac	hment A - Representative chromatograms	18

The chromatograms are graphical images are generated in the SCIEX MultiQuant software summary report. There is a chromatogram for each of the pentobarbital transitions. The first column is the pentobarbital quantitation ion transition (most abundant ion) chromatogram (m/z 225→182), identified as Pentobarbital 1. The second column is the chromatogram for the second most abundant ion transition (m/z 225→85), identified as Pentobarbital 2 and is used as a qualifier ion transition. The third column is the least abundant ion (m/z 225→ 138), identified as Pentobarbital 3 and is used as a qualifier transition. For the bone meal chromatograms (Figures A10 and A11, a fourth transition is included. It is directly below the Pentobarbital 1 figure. It is identified as Pentobarbital 4 and is the chromatogram for the transition m/z 225→ 42. It is more abundant than the pentobarbital quant transition and can be used as an alternative qualifier ion in cases of matrix interference peaks.

FOOD AND DRUG ADMINISTRATION OFFICE OF REGULATORY AFFAIRS Denver Laboratory	Document Number: WI-000140 Compendium method C-005.xx	Revision #: 04 Revised: Nov. 2019
Title: Determination of pentobarbital in ingredients of animal or foods using liquid chromatography tandem mass		Page 2 of 34

1. Purpose

This work instruction describes the application of LIB 4648 for the determination of pentobarbital in ingredients of animal origin and in finished pet foods using liquid chromatography tandem mass spectrometry.

2. Scope

For routine analysis of regulatory samples of numerous matrices of animal origin including: tallow (animal fat/grease), wet (canned) dog and cat food, dry dog and cat food, meat/bone meal, and meats including horse, beef, lamb, pork, goat, venison, bison, elk, rabbit, liver, and kidney using LC-MS/MS. Complete method performance details and validation data is summarized in the FDA QMiS Denver Lab (DENL) validation summary reports (RPRT- 00055, RPRT- 000060, RPRT- 00089, RPRT-000125, RPRT-000126).

3. References

Determination of Pentobarbital in Pet Food Using Liquid Chromatography Tandem Mass Spectrometry, Tuner, James A., Mohrhaus, Angie S., FCC SOP T064 Version 1. 3/1/2018.

- A. LC-HRMS Screen CALIFORNIA ANIMAL HEALTH &FOOD SAFETY LABORATORY, DTOX-02-877 Rev. 6. November 3, 2017.
- B. Pentobarbital in Beef Tallow March 2018, Personnel Correspondence with UC Davis, "Pentobarbital in Beef Tallow March 2018 (004) UC Davis extraction March 2018" I
- C. U.S. Food and Drug Administration, Office of Foods, Guidelines for the Validation of Chemical Methods for the FDA Foods Program, 2nd Edition 2015. https://www.fda.gov/downloads/scienceresearch/fieldscience/ucm27341
 - https://www.fda.gov/downloads/scienceresearch/fieldscience/ucm27341 8.pdf
- D. ORA Laboratory Manual, Volume II-Methods, Method Verification and Validation ORA-LAB.5.4.5 https://www.fda.gov/downloads/ScienceResearch/FieldScience/Laborat-oryManual/UCM092147.pdf
- E. DENL QMS # 18-3 and QMiS RPRT-000060 (Determination of pentobarbital in Tallow, wet pet food, and dry pet food using liquid chromatography tandem mass spectrometry).

FOOD AND DRUG ADMINISTRATION OFFICE OF REGULATORY AFFAIRS Denver Laboratory Title: Determination of pentobarbital in ingredients of animal origin and in finished pet foods using liquid chromatography tandem mass spectrometry. Document Number: WI-000140 Compendium method C-005.xx Revision #: 04 Revised: Nov. 2019 Page 3 of 34

- F. DENL QMS #18-4 and QMiS RPRT-000055 (determination of pentobarbital in tallow, wet pet food, and dry pet food using liquid chromatography tandem mass spectrometry).
- G. DENL QMS #18-4a and QMiS RPRT-000089 (determination of pentobarbital in wet dog food, and dry dog food using liquid chromatography tandem mass spectrometry).
- H. DENL QMS #18-4b and QMiS RPRT-000125(determination of pentobarbital in bone meal using liquid chromatography tandem mass spectrometry).
- DENL QMS #18-4c and QMiS RPRT-0000126 (determination of pentobarbital in meats, wet cat food, and dry cat food using liquid chromatography tandem mass spectrometry).
- J. U.S. Food and Drug Administration (2003) Guideline for Industry: Mass Spectrometry for Confirmation of the Identity of Animal Drug Residues, Fed. Regist. 68, 25617–25618. https://www.fda.gov/downloads/AnimalVeterinary/GuidanceComplianceEnforcement/GuidanceforIndustry/ucm052658.pdf
- K. CVM # GFI 118. Guidance for Industry Mass Spectrometry for Confirmation of the Identity of Animal Drug Residues https://www.fda.gov/media/70154/downloadhttp://www.fda.gov/downloads/ AnimalVeterinary/GuidanceComplianceEnforcement/GuidanceforIndustry/UCM0526 58.pdf
- L. DENL QMS #18-4a and QMiS RPRT-000089 (determination of pentobarbital in horsemeat, wet pet food, and dry pet food using liquid chromatography tandem mass spectrometry).\
- M. LIB 4648: Determination of pentobarbital in tallow using liquid chromatography tandem mass spectrometry (LC-MS/MS) CARTS: IR01702

http://inside.fda.gov:9003/downloads/PolicyProcedures/Laboratories/LaboratoryInformationBulletins/UCM632922.pdf

4. Procedure

4.1. Summary of Method

Pentobarbital is extracted from a homogenous portion of sample using acetonitrile. After centrifugation, the supernatant is diluted 1:1 with water and analyzed via LC-MS/MS using a multi-point solvent standard curve. Deuterated pentobarbital (pentobarbital-D5) is used as an internal standard

FOOD AND DRUG ADMINISTRATION OFFICE OF REGULATORY AFFAIRS Denver Laboratory	Document Number: WI-000140 Compendium method C-005.xx	Revision #: 04 Revised: Nov. 2019
Title: Determination of pentobarbital in ingredients of animal origin and in finished pet foods using liquid chromatography tandem mass spectrometry.		Page 4 of 34

(I.S.) to correct for sample matrix suppression and/or loss of analyte. Identification of pentobarbital in a sample is based on both correlation of pentobarbital chromatographic retention time (RT) with that of a standard(s) and ion ratio match.

4.2. Preparation of Samples

- 4.2.1. For tallow, at minimum a 25-gram sample portion is necessary. If sample appears heterogeneous, stir the 25 g sample manually with a spatula or spoon to ensure homogeneity.
- 4.2.2. For meat, at minimum a 50-gram sample portion is necessary. Tissue should be cut into small pieces, trying to avoid including any overly fatty portions, and then ground with dry ice in an appropriate blender. Resulting composite should be a fine powder.
- 4.2.3. For canned pet food, an entire can (approximately 150 grams or more) of the canned pet food is necessary. Wet pet food should be ground and mixed in in an appropriate blender. Do not add water. Dry ice may be used if necessary.
- 4.2.4. For dry pet food, approximately 150 grams is necessary. Dry pet food should be ground and mixed in an appropriate blender. Dry ice may be used if necessary (effective for products containing chewy 'bits').
- 4.2.5. For bone/meat meal, at minimum 50 grams is necessary. Bone meal should be ground and mixed in an appropriate blender. Dry ice may be used if necessary.
- 4.2.6. If dry ice is used in sample preparation, allow carbon dioxide gas to sublime in a freezer prior to removing aliquot for analysis.
- 4.2.7. If a sample appears heterogeneous, it should be mixed manually, or via a blender to ensure a homogeneous sample prior to removing a portion for analysis.
- 4.2.8. A 2.00-gram portion is used for each analysis. Two portions of each sample should be analyzed.

FOOD AND DRUG ADMINISTRATION OFFICE OF REGULATORY AFFAIRS Denver Laboratory	Document Number: WI-000140 Compendium method C-005.xx	Revision #: 04 Revised: Nov. 2019
Title: Determination of pentobarbital in ingredients of animal or foods using liquid chromatography tandem mass		Page 5 of 34

4.3. Preparation of Standards

- 4.3.1. Pentobarbital and pentobarbital-D5 are ordered premade with concentration 1000 μg/mL (1 mg/mL) in methanol.
- 4.3.2. Prepare a pentobarbital standard at 2,500 ng/mL with 25 μL of the 1000 μg/mL stock diluted to 10.0 mL with acetonitrile.
- 4.3.3. Prepare a pentobarbital ICV standard at 2,500 ng/mL with 25 μL of the 1000 μg/mL stock, from a different source or different ampoule, diluted to 10.0 mL with acetonitrile.
- 4.3.4. Prepare a pentobarbital-D5 internal standard spiking solution at 2,500 ng/mL with 25 μL of the 1000 μg/mL pentobarbital-D5 stock diluted to 10.0 mL with acetonitrile.
- 4.3.5. All prepared solutions have an expiration date of one year from the preparation date when stored at 4°C. If a component of the solution expires prior to one year of the preparation date, the earliest expiry date becomes the solution's expiration date.
- 4.3.6. Table 4.3.8 and 4.3.9 are examples of the preparations of working solutions and the solvent calibrants. Table 4.3.10 demonstrates the equivalent concentrations of the calibrants to the concentration in the samples for use in the processing method.
- 4.3.7. All eight calibration standards shown in Table 4.3.10 are not required for regulatory sample analysis; however, a minimum of 5 calibration standards are used with every batch of samples with a 10.0 ng/g standard (in sample amount), which is equivalent to an in vial concentration of 1.0 ng/mL as the lower limit of quantitation (LLOQ).

4.3.8. Working Solution Preparation (in acetonitrile)

Standard Name	starting concentration (μg/mL)	Standard added (mL)	Final vol (mL)	Final concentration (ng/mL)
Pentobarbital-2500	1,000	0.025	10.0	2,500
Pentobarbital ICV	1,000	0.025	10.0	2,500
Pent-D5 Spiking (ISTD)	1,000	0.025	10.0	2,500

FOOD AND DRUG ADMINISTRATION OFFICE OF REGULATORY AFFAIRS Denver Laboratory Title: Determination of pentobarbital in ingredients of animal origin and in finished pet foods using liquid chromatography tandem mass spectrometry. Document Number: WI-000140 Compendium method C-005.xx Revision #: 04 Revised: Nov. 2019 Page 6 of 34

4.3.9. Solvent Calibration Standard Preparation (in 50:50 acetonitrile:water)

Calibration	Initial conc	volume of	Volume D5-	Final	Final Conc	Final Conc
Curve	pentobarbital	pentobarbital std	ISTD Added	Volume	pentobarbital	d5-Pent
Curve	(ng/mL)	added (mL)	(2500 ng/mL)	(mL)	(ng/mL)	(ng/mL)
Cal-1	2,500	0.020	0.100	50.0	1.00	5.00
Cal-2	2,500	0.010	0.040	20.0	1.25	5.00
Cal-3	2,500	0.010	0.020	10.0	2.50	5.00
Cal-4	2,500	0.020	0.020	10.0	5.00	5.00
Cal-5	2,500	0.040	0.020	10.0	10.00	5.00
Cal-6	2,500	0.100	0.020	10.0	25.00	5.00
Cal-7	2,500	0.200	0.020	10.0	50.00	5.00
Cal-8	2,500	0.400	0.020	10.0	100.00	5.00
ICV	2,500	0.040	0.020	10.0	10.00	5.00

4.3.10. Conversion of Solvent Standards to In Sample Concentration

Calibration Curve	In Vial Final Concentratio n (ng/mL)	Sample weight. (g)	Volume of ACN extraction (mL)	Dilution	Equivalent pentobarbital in matrix concentration (ng/g)	Equivalent d5-pent (ISTD) In matrix concentration (ng/g)
Cal-1	1.00	2.00	10	2.00	10.0	50.0
Cal-2	1.25	2.00	10	2.00	12.5	50.0
Cal-3	2.50	2.00	10	2.00	25.0	50.0
Cal-4	5.00	2.00	10	2.00	50.0	50.0
Cal-5	10.0	2.00	10	2.00	100	50.0
Cal-6	25.0	2.00	10	2.00	250	50.0
Cal-7	50.0	2.00	10	2.00	500	50.0
Cal-8	100.0	2.00	10	2.00	1000	50.0
ICV	10.0	2.00	10	2.00	100	50.0

FOOD AND DRUG ADMINISTRATION OFFICE OF REGULATORY AFFAIRS Denver Laboratory	Document Number: WI-000140 Compendium method C-005.xx	Revision #: 04 Revised: Nov. 2019
Title: Determination of pentobarbital in ingredients of animal or foods using liquid chromatography tandem mass		Page 7 of 34

4.3.11. Fortified (spiked) sample preparation (demonstrates typical fortification levels for pentobarbital in sample)

Spike Level (ng/g)	initial concentration (ng/mL)	volume used (mL)	sample weight (g)	final concentration in sample (ng/g)
10 ng/g	2500	0.008	2.00	10.0
50 ng/g	2500	0.040	2.00	50.0
250 ng/g	2500	0.200	2.00	250

4.3.1. Example spiking calculation:

40 μL	2500 ng	= 50.0 ng
2.00 grams of negative control	mL	g

4.3.2. Example calculation of in vial versus in sample concentration

50 ng	2 g sample amount	0.500 mL extract	= 5.00 ng in vial
g	10 mL extraction volume	1.00 mL final volume	mL

FOOD AND DRUG ADMINISTRATION OFFICE OF REGULATORY AFFAIRS Denver Laboratory	Document Number: WI-000140 Compendium method C-005.xx	Revision #: 04 Revised: Nov. 2019
Title: Determination of pentobarbital in ingredients of animal or foods using liquid chromatography tandem mass		Page 8 of 34

4.4. Reagents

- 4.4.1. Water, Fisher, LC-MS grade
- 4.4.2. Acetonitrile, Fisher, LC-MS Grade
- 4.4.3. Pentobarbital, Cerilliant, 1.000 ± 0.005 mg/mL in methanol, 1 mL ampoule, part # P-010
- 4.4.4. Pentobarbital-D₅, Cerilliant, 1.000 ± 0.005 mg/mL in methanol, 1 mL ampoule, part #: P-013
- 4.4.5. Diluent for standards, 50/50 water/acetonitrile (v/v), 500 mL. Combine 250 mL water + 250 mL acetonitrile in a 500 mL graduated cylinder. Cap and invert to mix.
- 4.5. Equipment (equivalent equipment may be substituted)
 - 4.5.1. Vortexer/ Mixer, Troemner, (500-2500 rpm)
 - 4.5.2. SPEX Geno/Grinder 2000 (500 rpm)
 - 4.5.3. Sonicator (Branson 2510 or 8510)
 - 4.5.4. Appropriate mixers, blenders, food processors, etc. used to homogenize sample matrix if necessary
 - 4.5.5. Centrifuge capable of 6000 rpm with refrigeration (4 °C) for 50 mL tubes
 - 4.5.6. Plastic centrifuge tubes with caps, 15 mL and 50 mL
 - 4.5.7. Microcentrifuge tubes, at least 1 mL capacity
 - 4.5.8. Nylon syringe filters, PALL Life Science Acrodisc 13 mm Syringe Filters 0.2 μm
 - 4.5.9. Luer slip 1 mL syringes
 - 4.5.10. 2 mL glass amber autosampler vials and pre-slit snap caps (#66030-608)
 - 4.5.11. Calibrated pipettes and volumetric glassware

FOOD AND DRUG ADMINISTRATION OFFICE OF REGULATORY AFFAIRS Denver Laboratory	Document Number: WI-000140 Compendium method C-005.xx	Revision #: 04 Revised: Nov. 2019
Title: Determination of pentobarbital in ingredients of animal of foods using liquid chromatography tandem mass		Page 9 of 34

- 4.6. Instrumentation (equivalent instrumentation may be substituted)
 - 4.6.1. Agilent 1200 HPLC system, Combi Pal autosampler, with an AB SCIEX 5500 QTRAP MS instrument operated in the negative mode with an ESI source using Multiple Reaction Monitoring (MRM). AB SCIEX 1.6.2 software was used for instrument control and MultiQuant 3.0 was used for data processing.
 - 4.6.2. AB SCIEX QTRAP 5500 source/parameter settings:

A. Curtain gas: 30 psi

B. GS1: 50 psiC. GS2: 60 psi

D. Collision gas: medium

E. Ion spray voltage (IS): -3500VF. Source temperature: 400°CG. Entrance Potential (EP): -10V

4.6.3. Pentobarbital MS parameters: Retention times (RT), transitions, declustering potential (DP), collision energy (CE), cell exit potential (CXP), and the resulting typical ion ratios for the product ions of each analyte from the ABI SCIEX 5500 QTRAP analysis.

	Typical RT (min)	1	ransitic	on	ISTD	DP (V)	CE (V)	CXP (V)	Average ion ratio, qual/quant %
Analyte	()		(/			(-)	(-)	(-)	
				182			-19	-13	100
pentobarbital	4.20	225	$\bar{5} \longrightarrow$	85	Pent-D5	-100	-18	-9	15
periobarbitai	7.20	220		138	T CIR DO		-21	-10	7
				42*			-46	-7	275*
pentobarbital-D5	4.20	230	\rightarrow	187		-100	-17	-10	N/A

^{*}additional transition of 225→42 was added as additional/alternative confirmation criteria due to matrix interference peak for 225→85 transition in some bone meal sources at <10 ppb level. The 225→42 transition is more abundant than the 225→82 transition, giving the ion ratio of >100%.

FOOD AND DRUG ADMINISTRATION OFFICE OF REGULATORY AFFAIRS Denver Laboratory	Document Number: WI-000140 Compendium method C-005.xx	Revision #: 04 Revised: Nov. 2019
Title: Determination of pentobarbital in ingredients of animal o foods using liquid chromatography tandem mass		Page 10 of 34

4.6.4. Combi Pal autosampler settings

A. sample injection volume: 5 μL

B. autosampler tray temperature: 15°C

C. Combi Pal Injector wash solution 1: 95% water/5% acetonitrile

D. Combi Pal Injector wash solution 2: 95% acetonitrile/5% water

4.6.5. Agilent 1200 HPLC settings

- A. HPLC column Agilent Zorbax Eclipse Plus C18, 2.1 x 50 mm, 1.8 micron size column (part # 959757-902)
- B. A divert valve directed column effluent to waste before (0-0.5 minutes), and during the system re-equilibration time (5.0-8.50)
- C. Column temperature was 40°C
- D. LC flow rate was 0.350 mL/min.
- E. The mobile phase was water (A) and acetonitrile (B), and the LC gradient is described in Table 4.6.5.F below
- F. LC gradient for pentobarbital

@Step	Total Time (min)	Flow Rate (µl/min)	A (%) (Water)	B (%) (Acetonitrile)
0	0.00	350	95.0	5.0
1	3.50	350	5.0	95.0
2	4.50	350	95.0	5.0
4	8.50	350	95.0	5.0

LC note: It has been observed that any mobile phase modifiers (formic acid, ammonium formate, etc) have a significant negative effect on pentobarbital response. Ensure LC lines have been thoroughly cleaned and purged with water and acetonitrile before analysis.

FOOD AND DRUG ADMINISTRATION OFFICE OF REGULATORY AFFAIRS Denver Laboratory Title: Determination of pentobarbital in ingredients of animal origin and in finished pet foods using liquid chromatography tandem mass spectrometry. Pocument Number: WI-000140 Compendium method C-005.xx Revision #: 04 Revised: Nov. 2019 Page 11 of 34

4.7. Extraction

- A. Weigh 2.00± 0.05 g of each homogenized sample into a 50 mL centrifuge tube. For each unknown sample, weigh out two portions.
- B. For each batch, include an empty tube to serve as Reagent Blank (RB).
- C. Weigh out three portions of negative control material to serve as negative control (NC), matrix spike (SPK), and matrix spike duplicate (DUP).
- D. For all samples in the batch, including RB, NC, SPK, and DUP, add 40 μ L of 2500 ng/mL pentobarbital-D5, resulting in a 50 ng/g d5-pentobarbital concentration.
- E. Fortify spike (SPK) and duplicate (DUP) portions with 40 μL of 2500 ng/mL pentobarbital spiking standard, resulting in a 50 ng/g pentobarbital concentration in sample. (Fortification level may be adjusted as necessary, as long as the concentration falls within the calibration curve).
- F. If a sample appears to be significantly different than an available control matrix, an additional portion of sample should be spiked to assess extraction efficiency and matrix effects compared to the most similar control matrix.
- G. Add 10 mL of acetonitrile to each sample tube.
- H. Cap tubes tightly and shake @ 500 rpm for 5 minutes on the geno grinder.
- I. Sonicate sample tubes for 30 minutes.
- J. Vortex sample tubes for 30 seconds.
- K. Centrifuge sample tubes at 6000 x g and 4°C for 10 minutes.
- L. Combine 500μL of sample supernatant with 500μL of water in a microcentrifuge tube; vortex to mix, then filter using a 0.2 μm Nylon syringe filter into a 2 mL amber LC vial and cap.
- M. Analyze via LC-MS/MS.

FOOD AND DRUG ADMINISTRATION OFFICE OF REGULATORY AFFAIRS Denver Laboratory	Document Number: WI-000140 Compendium method C-005.xx	Revision #: 04 Revised: Nov. 2019
Title: Determination of pentobarbital in ingredients of animal or foods using liquid chromatography tandem mass		Page 12 of 34

4.8. Data Analysis and Quality Acceptance

- 4.8.1. Calibration curves are established from a multi-point solvent calibrant standard curve, typically ranging from 1-50 ng/mL in vial, equivalent 10-500 ng/g in sample, with the concentration on the x-axis and internal standard corrected peak response on the y-axis. Suggested seven-point curve for routine sample analysis is 1.00, 1.25, 2.5, 5.0, 10.0, 25.0, 50.0 ng/mL (equivalent to 10.0, 12.5, 25, 50, 100, 250, 500 ng/g in sample).
- 4.8.2. Quantitative results will be reported for samples with responses that fall within the standard curve range and meet identity confirmation criteria.
- 4.8.3. The calculated method limit of detection (MDL) and calculated method limit of quantitation (LOQ), and typical linearity (r²) are indicated in Table 4.8.4.

FOOD AND DRUG ADMINISTRATION OFFICE OF REGULATORY AFFAIRS Denver Laboratory	Document Number: WI-000140 Compendium method C-005.xx	Revision #: 04 Revised: Nov. 2019
Title: Determination of pentobarbital in ingredients of animal or foods using liquid chromatography tandem mass		Page 13 of 34

4.8.4. MDL, LOQ, Linearity Table

Matrix	MDL (ng/g) n=7*	LOQ (ng/g) n=7*	Average Linearity (r²)
	Finished Pet F	ood	
Dry Dog Food	2.9	9.3	0.9989
Wet Dog Food	2.5	7.8	0.9994
Dry Cat Food	1.8	5.8	0.9999
Wet Cat Food	1.0	3.2	1.0000
	Rendering Stre	ams	
Tallow	3.6	11.6	0.9991
Bone Meal * n=9	2.3	8.0	0.9991
	Meat Produc	ts	
Beef	1.6	5.2	0.9997
Bison	1.0	3.2	0.9998
Elk	1.2	3.9	0.9994
Goat	1.3	4.2	0.9998
Horse	0.9	2.9	0.9991
Lamb	0.9	3.0	0.9998
Pork	0.9	3.0	1.0000
Rabbit	1.6	5.2	0.9999
Venison	0.8	2.6	1.0000
Kidney (Bison)	1.8	5.7	0.9999
Liver (Beef)	1.6	5.0	0.9997
Average for all matrices:	1.6	5.3	0.9996

FOOD AND DRUG ADMINISTRATION OFFICE OF REGULATORY AFFAIRS Denver Laboratory	Document Number: WI-000140 Compendium method C-005.xx	Revision #: 04 Revised: Nov. 2019
Title: Determination of pentobarbital in ingredients of animal or foods using liquid chromatography tandem mass		Page 14 of 34

- 4.8.5. The tallow MDL and LOQ data in LIB 4648 were originally calculated from 12.5 ng/g spikes (n=9), resulting in calculated MDL of 2.4 ng/g and LOQ of 8.2 ng/g). Additional tallow spikes (n=7) were analyzed at 5 ng/g and 10 ng/g and the MDL and LOQ were re-calculated using that data, which is in Table 4.8.4 above.
- 4.8.6. All calibration curves were generated with the AB SCIEX MultiQuant software. Depending on the range of the calibration curve, a linear fit with no weighting or with 1/x weighting (not forced through zero) can be used for all recovery calculations. If a smaller dynamic range is used, a linear curve with no weighting should be used. All calibration curves should have r² ≥0.995.
- 4.8.7. Precision and accuracy general guidance (performance may vary with sample matrix, especially for different tallow/animal fat matrices):
 - A. The FDA OFVM specifies that analyte recovery should be within the range 60%-115% corresponding to concentration from 10-100 mg/kg (ppb), with RSD_r \leq 22%. Recoveries for elk (123%) and beef liver (121%) were >115% at the 10 ng/g spiking level. Higher recoveries at the low spiking level may be due to the use of a linear curve without weighting.
 - B. The aforementioned ranges are for guidance on data acceptability. Ranges for precision and accuracy should be determined in house after a sufficient number of data points are obtained.
 - C. Reagent blank (RB) and negative control (NC) should have responses below the lowest calibration standard.
 - D. Validation data demonstrated satisfactory quantitative analysis for pentobarbital determined in all matrices at spiking levels 10, 50, and 250 ng/g, with method accuracy generally ranging from 85-123%, with RSD_r \leq 16%.
 - E. Any QC failures must be investigated prior to reporting results.

FOOD AND DRUG ADMINISTRATION OFFICE OF REGULATORY AFFAIRS Denver Laboratory Title: Determination of pentobarbital in ingredients of animal origin and in finished pet foods using liquid chromatography tandem mass spectrometry. Document Number: WI-000140 Compendium method C-005.xx Revision #: 04 Revised: Nov. 2019 Page 15 of 34

- F. If a sample demonstrates a response above the highest calibration point the sample can be further diluted with the 50:50 acetonitrile:water diluent to approximately the midpoint of the solvent curve. The vials (calibration curve, quality control samples, and diluted sample(s)) can then be reinjected and the data can be processed without the internal standard correction. Appropriateness of this approach should be evaluated on a case by case basis. Care must be taken to ensure the concentration range of the solvent curve reflects the dilution of the sample. (Data was evaluated with and without the use of internal standard during validation for some matrices). It is advised to use the internal standard correction for extracts that fall within the routine calibration curve range.
- 4.8.8. Positive confirmation of identity for positive samples, and spiked samples
 - A. Signal to noise must be >3:1 (The AB SCIEX MultiQuant software is used to calculate signal to noise, if required.)
 - B. Retention time must match the comparison standard(s) within 5%.
 - C. Ion ratios must match the comparison standard(s) by an absolute value of ±20%. (MultiQuant software uses the average ion ratio of all standards in the calibration curve). If the ratio of either of the qualifier transitions fails due to instrument sensitivity, the other two transitions may be used with a ± 10% tolerance (absolute) on the ratio when compared to the average standard ratio.

4.8.9. Sample Result Reporting

- A. Quantitative sample results greater than the lower limit of quantitation (LLOQ) of 10 ng/g shall be reported.
- B. Samples with calculated amounts less than the lower limit of quantitation (LLOQ) of 10 ng/g, but greater than the MDL shall be reported as "pentobarbital detected at <10 ng/g, but greater than MDL", along with the value for the MDL for that matrix.
- C. Samples with calculated amounts < MDL shall be reported as "pentobarbital not detected at or above MDL".

FOOD AND DRUG ADMINISTRATION OFFICE OF REGULATORY AFFAIRS Denver Laboratory Title: Determination of pentobarbital in ingredients of animal origin and in finished pet foods using liquid chromatography tandem mass spectrometry. Pocument Number: Revision #: 04 Revised: Nov. 2019 Page 16 of 34

5. Glossary

- A. RB: Reagent Blank. Used to verify reagents are uncontaminated by interfering components, the reagent blank is an extract that contains no sample matrix. Carried thorough the extraction as if it were a sample, one must be extracted with each batch and display no interference peaks at the reference times of interest at or above lowest calibration.
- H. NC: Negative Control. Used to verify the lack of matrix effects, the control is an aliquot of matrix material known to contain no analytes of interest. One must be extracted with each batch for each matrix type, and must display no interference peaks at reference times of interest at or above lowest calibration.
- I. SPK/DUP: Matrix spike/matrix spike duplicate. Used to demonstrate effective and reproducible extraction, the matrix spike and duplicate are two aliquots of negative control matrix material, each fortified at a level near the midpoint of the curve. A pair of matrix spikes must be extracted and analyzed with each batch for each matrix type.
- J. ICV: Independent Calibration Verification. Used to assure the accuracy of the calibration curve, the ICV is a solvent standard prepared from a secondary standard source.
- K. CCV: Continuing Calibration Verification. Used to check the calibration during a run, the CCV is a re-injection of a midpoint solvent standard curve. A CCV is injected after every ten sample extracts and at the end of the analytical sequence.
- L. MDL = Method Detection Limit = $\sigma(t_{(df=N-1,1-\alpha=0.99,one\ sided)})$
- M. Where: σ = standard deviation (of 5 ng/g spikes, N=7)
- N. t = Student's T value for df = N-1 at the 99% confidence level, one sided
- O. LOQ = Limit of Quantitation = $\sigma(10)$
- P. Where: σ = standard deviation (of 10 ng/g spikes, N=7)
- Q. LLOQ = Lower Limit of Quantitation, lowest standard in calibration curve. For this analysis, the LLOQ should be equal to 10 ng/g in sample (or lower).

6. Records

- A. Analyst Worksheets
- B. Batch Records

FOOD AND DRUG ADMINISTRATION OFFICE OF REGULATORY AFFAIRS Denver Laboratory	Document Number: WI-000140 Compendium method C-005.xx	Revision #: 04 Revised: Nov. 2019
Title: Determination of pentobarbital in ingredients of animal or foods using liquid chromatography tandem mass		Page 17 of 34

C. Equipment FVPMS

7. Document History

Revision #	Status* (D, I, R)	Date	Author Name and Title	Approving Official Name and Title
00	I	11/30/2018	Tara Nickel, Chemist Christine Casey, Chemist	R. Stadtmuller, Quality System Manager
01	R	12/4/18	Tara Nickel, Chemist	R. Stadtmuller, Quality System Manager
02	R	03/06/19	Tara Nickel, Chemist	R. Stadtmuller, Quality System Manager
03	R	5/1/2019	Tara Nickel, Chemist	R. Stadtmuller, Quality System Manager
04	R	11/13/2019	T. Nickel, Chemist	R. Stadtmuller, Quality System Manager

^{* -} D: Draft, I: Initial, R: Revision

8. Change History

Revision #	Change
00	Original
01	Change 10 uL to 20 uL on 2.7.2. Add reporting statement on 2.8.9.3 for amounts below 10ng/g and >MDL.
02	Minor corrections for spelling, grammar, and formatting
03	Add dry dog food, wet dog food, horsemeat information throughout document
04	Add information for additional matrices throughout document. Copy to ORS template.

9. Attachments

Attachment A - Representative chromatograms	. 18-29
Attachment B – Instrument Acquisition method	. 30-34

FOOD AND DRUG ADMINISTRATION OFFICE OF REGULATORY AFFAIRS Denver Laboratory	Document Number: WI-000140 Compendium method C-005.xx	Revision #: 04 Revised: Nov. 2019
Title: Determination of pentobarbital in ingredients of animal o foods using liquid chromatography tandem mass	Page 18 of 34	

Attachment A - Representative chromatograms


The result summary table states the analyte peak precursor name, observed analyte retention time (RT), expected analyte retention time (RT) time, calculated concentration (ng/g), analyte response (peak area), calculated Ion Ratio (ration of qualifier ion peak area to quantifier ion peak area), and Ratio Confirmation (a checkmark indicates the ratio matches the comparison standards ±20%).

The chromatograms are graphical images are generated in the SCIEX MultiQuant software summary report. There is a chromatogram for each of the pentobarbital transitions. The first column is the pentobarbital quantitation ion transition (most abundant ion) chromatogram (m/z $225\rightarrow182$), identified as Pentobarbital 1. The second column is the chromatogram for the second most abundant ion transition (m/z $225\rightarrow85$), identified as Pentobarbital 2 and is used as a qualifier ion transition. The third column is the least abundant ion (m/z $225\rightarrow138$), identified as Pentobarbital 3 and is used as a qualifier transition. For the bone meal chromatograms (Figures A10 and A11, a fourth transition is included. It is directly below the Pentobarbital 1 figure. It is identified as Pentobarbital 4 and is the chromatogram for the transition m/z $225\rightarrow42$. It is more abundant than the pentobarbital quant transition and can be used as an alternative qualifier ion in cases of matrix interference peaks.

FOOD AND DRUG ADMINISTRATION OFFICE OF REGULATORY AFFAIRS Denver Laboratory	Document Number: WI-000140 Compendium method C-005.xx	Revision #: 04 Revised: Nov. 2019
Title: Determination of pentobarbital in ingredients of animal or foods using liquid chromatography tandem mass	Page 19 of 34	

Figure A1 Result Summary for: Solvent Calibrant # 3- equivalent to ng/g in sample (In vial concentration: 5.0 ng/mL pentobarbital, 5.0 ng/mL d5-pentobarbital)


Analyte Peak Name	Analyte RT	Expected RT	Calculated Concentration (ng/g)	Analyte Response	Calculated Ion Ratio (Expected Value)	Ratio Confirms
Pentobarbital 1 (225->182.0)	4.21	4.20	52.50	170000.0		
Pentobarbital 2 (225->85.0)	4.21	4.20		26600	15.6% (15.8%)	√
Pentobarbital 3 (225->138.0)	4.21	4.20		11500	6.8% (7.2%)	√

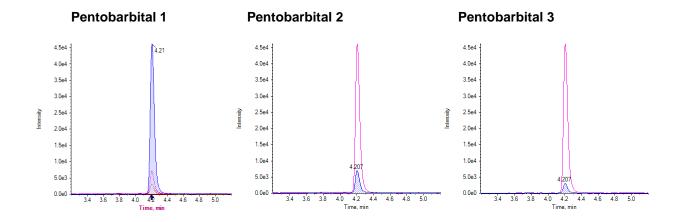

FOOD AND DRUG ADMINISTRATION OFFICE OF REGULATORY AFFAIRS Denver Laboratory	Document Number: WI-000140 Compendium method C-005.xx	Revision #: 04 Revised: Nov. 2019
Title: Determination of pentobarbital in ingredients of animal or foods using liquid chromatography tandem mass		Page 20 of 34

Figure A2: Result Summary for Tallow Negative Control, Source T1


Analyte Peak Name	Analyte RT	Expected RT	Calculated Concentration (ng/g)	Analyte Response	Calculated Ion Ratio (Expected Value)	Ratio Confirms
Pentobarbital 1 (225->182.0)	3.90	4.20	< 0	3190.0		
Pentobarbital 2 (225->85.0)	4.22	4.20		1920	60.1% (15.8%)	
Pentobarbital 3 (225->138.0)	4.06	4.20		395	12.4% (7.2%)	√

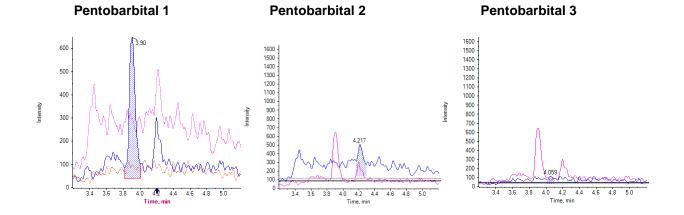

FOOD AND DRUG ADMINISTRATION OFFICE OF REGULATORY AFFAIRS Denver Laboratory	Document Number: WI-000140 Compendium method C-005.xx	Revision #: 04 Revised: Nov. 2019
Title: Determination of pentobarbital in ingredients of animal or foods using liquid chromatography tandem mass	Page 21 of 34	

Figure A3: Result Summary for Tallow - Fortified at 50 ng/g


Analyte Peak Name	Analyte RT	Expected RT	Calculated Concentration (ng/g)	Analyte Response	Calculated Ion Ratio (Expected Value)	Ratio Confirms
Pentobarbital 1 (225_>182.0)	4.21	4.20	51.65	129000.0		
Pentobarbital 2 (225->85.0)	4.21	4.20		20500	15.8% (15.8%)	✓
Pentobarbital 3 (225->138.0)	4.21	4.20		8930	6.9% (7.2%)	✓

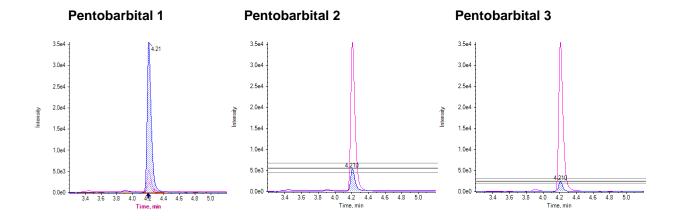

FOOD AND DRUG ADMINISTRATION OFFICE OF REGULATORY AFFAIRS Denver Laboratory	Document Number: WI-000140 Compendium method C-005.xx	Revision #: 04 Revised: Nov. 2019
Title: Determination of pentobarbital in ingredients of animal or foods using liquid chromatography tandem mass	Page 22 of 34	

Figure A4: Result Summary for Wet Dog Food Negative Control, Source W2


Analyte Peak Name	Analyte RT	Expected RT	Calculated Concentration (ng/g)	Analyte Response	Calculated Ion Ratio (Expected Value)	Ratio Confirms
Pentobarbital 1 (225->182.0)	3.90	4.20	<0	1326.0		
Pentobarbital 2 (225->85.0)	3.85	4.20		824	62.2% (14.3%)	
Pentobarbital 3 (225->138.0)	4.29	4.20		243	18.4% (6.8%)	✓

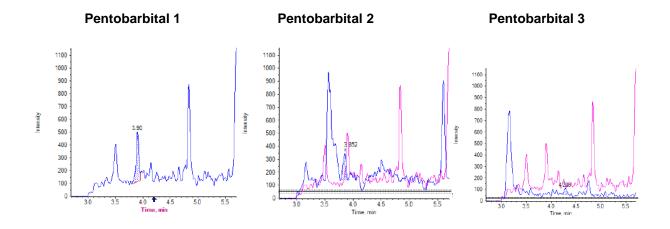

FOOD AND DRUG ADMINISTRATION OFFICE OF REGULATORY AFFAIRS Denver Laboratory	Document Number: WI-000140 Compendium method C-005.xx	Revision #: 04 Revised: Nov. 2019
Title: Determination of pentobarbital in ingredients of animal or foods using liquid chromatography tandem mass	Page 23 of 34	

Figure A5: Result Summary for Wet Dog Food Fortified at 50 ng/g, Source W2


Analyte Peak Name	Analyte RT	Expected RT	Calculated Concentration (ng/g)	Analyte Response	Calculated Ion Ratio (Expected Value)	Ratio Confirms
Pentobarbital 1 (225->182.0)	4.19	4.20	56.02	53007.0		
Pentobarbital 2 (225->85.0)	4.19	4.20		6812	12.9% (14.3%)	✓
Pentobarbital 3 (225->138.0)	4.19	4.20		3619	6.8% (6.8%)	✓

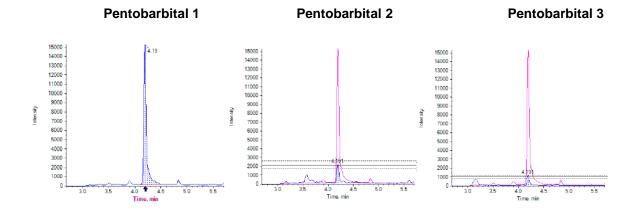

FOOD AND DRUG ADMINISTRATION OFFICE OF REGULATORY AFFAIRS Denver Laboratory	Document Number: WI-000140 Compendium method C-005.xx	Revision #: 04 Revised: Nov. 2019
Title: Determination of pentobarbital in ingredients of animal or foods using liquid chromatography tandem mass	Page 24 of 34	

Figure A6: Result Summary for Dry Dog Food Negative Control, Source D3


Analyte Peak Name	Analyte RT	Expected RT	Calculated Concentration (ng/g)	Analyte Response	Calculated Ion Ratio (Expected Value)	Ratio Confirms
Pentobarbital 1 (225->182.0)	N/A	4.20	N/A	1326.0	N/A	
Pentobarbital 2 (225->85.0)	3.95	4.20		824	0.0% (14.3%)	
Pentobarbital 3 (225->138.0)	3.92	4.20		243	0.0% (6.8%)	

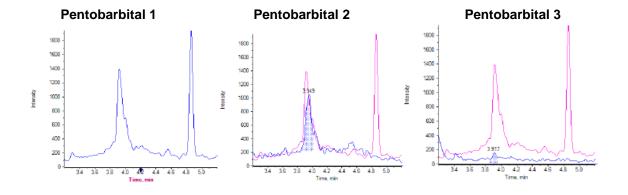

FOOD AND DRUG ADMINISTRATION OFFICE OF REGULATORY AFFAIRS Denver Laboratory	Document Number: WI-000140 Compendium method C-005.xx	Revision #: 04 Revised: Nov. 2019
Title: Determination of pentobarbital in ingredients of animal or foods using liquid chromatography tandem mass		Page 25 of 34

Figure A7: Result Summary for Dry Dog Food Fortified at 50 ng/g, Source D3


Analyte Peak Name	Analyte RT	Expected RT	Calculated Concentration (ng/g)	Analyte Response	Calculated Ion Ratio (Expected Value)	Ratio Confirms
Pentobarbital 1 (225->182.0)	4.22	4.20	51.26	68852.0		
Pentobarbital 2 (225->85.0)	4.22	4.20		9158	13.3% (14.3%)	✓
Pentobarbital 3 (225->138.0)	4.22	4.20		5022	7.3% (6.9%)	✓

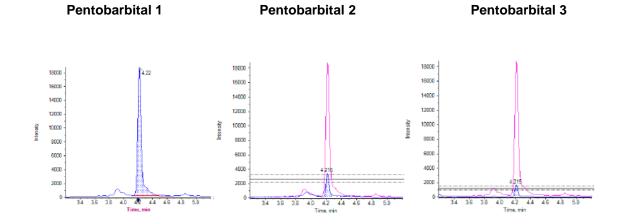

FOOD AND DRUG ADMINISTRATION OFFICE OF REGULATORY AFFAIRS Denver Laboratory	Document Number: WI-000140 Compendium method C-005.xx	Revision #: 04 Revised: Nov. 2019
Title: Determination of pentobarbital in ingredients of animal or foods using liquid chromatography tandem mass		Page 26 of 34

Figure A8: Result Summary for Horsemeat Negative Control


Analyte Peak Name	Analyte RT	Expected RT	Calculated Concentration (ng/g)	Analyte Response	Calculated Ion Ratio (Expected Value)	Ratio Confirms
Pentobarbital 1 (225->182.0)	N/A	4.69	N/A	N/A		
Pentobarbital 2 (225->85.0)	4.53	4.69		2165	0.0% (14.9%)	
Pentobarbital 3 (225->138.0)	4.54	4.69		780	0.0% (6.9%)	

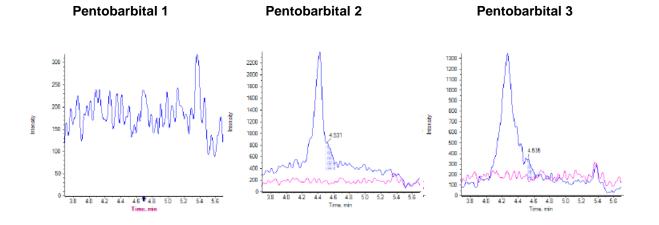

FOOD AND DRUG ADMINISTRATION OFFICE OF REGULATORY AFFAIRS Denver Laboratory	Document Number: WI-000140 Compendium method C-005.xx	Revision #: 04 Revised: Nov. 2019
Title: Determination of pentobarbital in ingredients of animal of foods using liquid chromatography tandem mas	Page 27 of 34	

Figure A9: Result Summary for Horsemeat Fortified at 50 ng/g

Analyte Peak Name	Analyte RT	Expected RT	Calculated Concentration (ng/g)	Analyte Response	Calculated Ion Ratio (Expected Value)	Ratio Confirms
Pentobarbital 1 (225->182.0)	4.69	4.69	50.29	70189.0		
Pentobarbital 2 (225->85.0)	4.53	4.69		11187	15.9% (14.9%)	✓
Pentobarbital 3 (225->138.0)	4.54	4.69		4374	6.2% (6.9%)	✓

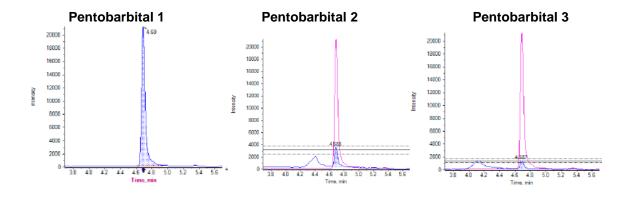


FOOD AND DRUG ADMINISTRATION OFFICE OF REGULATORY AFFAIRS Denver Laboratory	Document Number: WI-000140 Compendium method C-005.xx	Revision #: 04 Revised: Nov. 2019
Title: Determination of pentobarbital in ingredients of animal o foods using liquid chromatography tandem mass	Page 28 of 34	


Figure A10: Result Summary for Bone Meal Negative Control (source 2)

Analyte Peak Name	Analyte RT	Expecte d RT	Calc Conc. (ng/g)	Analyte Response	Calculated Ion Ratio (Expected Value)	Ratio Confirms
Pentobarbital 1 (225->182.0)	N/A	4.23	N/A	N/A		
Pentobarbital 2 (225->85.0)	4.25	4.23		7.07e3	0.0% (15.9%)	
Pentobarbital 3 (225->138.0)	N/A	4.23		N/A	0.0% (6.1%)	
Pentobarbital 4 (225->42.0)	4.23	4.23		5.24e3	0.0% (281.7%)	

Chromatograms – Bars on peaks are expected ion ratio ± 20% of comparison standards

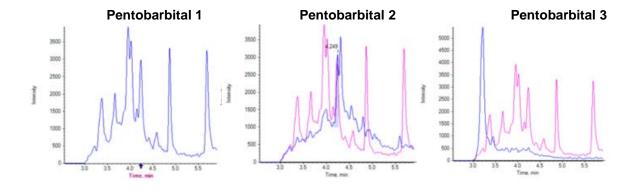

Pentobarbital 4

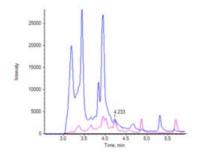

FOOD AND DRUG ADMINISTRATION OFFICE OF REGULATORY AFFAIRS Denver Laboratory	Document Number: WI-000140 Compendium method C-005.xx	Revision #: 04 Revised: Nov. 2019
Title: Determination of pentobarbital in ingredients of animal or foods using liquid chromatography tandem mass		Page 29 of 34


Figure A11: Result Summary for Bone Meal Fortified at 50 ng/g

Analyte Peak Name	Analyte RT	Expecte d RT	Calc Conc. (ng/g)	Analyte Response	Calculated Ion Ratio (Expected Value)	Ratio Confirms
Pentobarbital 1 (225->182.0)	4.25	4.23	45.76	1.93e5		
Pentobarbital 2 (225->85.0)	4.25	4.23		4.87e4	25.3% (15.9%)	√
Pentobarbital 3 (225->138.0)	4.24	4.23		1.54e4	8.0% (6.1%)	✓
Pentobarbital 4 (225->42.0)	4.24	4.23		5.16e5	267.4% (281.7%)	✓

Pentobarbital 2


Pentobarbital 3

Chromatograms – Bars on peaks are expected ion ratio ± 20% of comparison standards

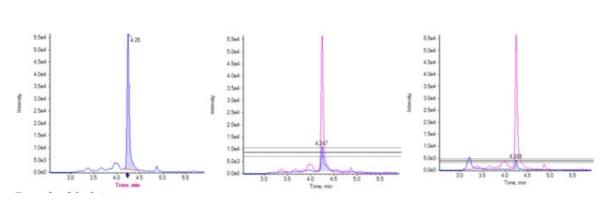

Pentobarbital 4

Pentobarbital 1

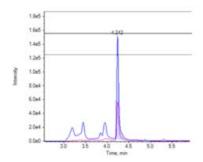

FOOD AND DRUG ADMINISTRATION OFFICE OF REGULATORY AFFAIRS Denver Laboratory Title: Determination of pentobarbital in ingredients of animal origin and in finished pet foods using liquid chromatography tandem mass spectrometry. Document Number: WI-000140 Compendium method C-005.xx Revision #: 04 Revised: Nov. 2019 Page 30 of 34

Attachment B: AB SCIEX QTRAP 5500 data acquisition method

HPLC column – Agilent Zorbax Eclipse Plus C18, 2.1 x 50 mm, 1.8 micron size column

Mobile Phase. A: Water

Mobile Phase. B: Acetonitrile

Synchronization Mode: LC Sync

Auto-Equilibration: Off

Acquisition Duration: 8min31sec

Number Of Scans: 1854

Periods In File: 1

Acquisition Module: Acquisition Method

Software version Analyst 1.6.2

MS Method Properties:

Period 1:

Scans in Period: 1854

Relative Start Time: 1000.00 msec

Experiments in Period: 1

Period 1 Experiment 1:

Scan Type: MRM (MRM)
Scheduled MRM: No
Polarity: Negative

Scan Mode: N/A

Ion Source: Turbo Spray

Resolution Q1: Unit
Resolution Q3: Unit
Intensity Thres.: 0.00 cps
Settling Time: 0.0000 msec
MR Pause: 5.0070 msec

MCA: No

FOOD AND DRUG ADMINISTRATION OFFICE OF REGULATORY AFFAIRS Denver Laboratory	Document Number: WI-000140 Compendium method C-005.xx	Revision #: 04 Revised: Nov. 2019
Title: Determination of pentobarbital in ingredients of animal or foods using liquid chromatography tandem mass	Page 31 of 34	

Step Size: 0.00 Da

@Q1 Mass (Da)Q3 Mass (Da) Dwell(msec) Param Start Stop ID

225.000 182.000 50.00 CE -19.00 -19.00 Pentobarbital 1

CXP -13.00 -13.00

@Q1 Mass (Da) Q3 Mass (Da) Dwell(msec) Param Start Stop ID 225.000 85.000 50.00 CE -18.00 -18.00 Pentobarbital 2 CXP -9.00-9.00

@Q1 Mass (Da) Q3 Mass (Da) Dwell(msec) Param Start Stop ID 225.000 138.000 50.00 CE -21.00 -21.00 Pentobarbital 3 CXP -10.00 -10.00

@Q1 Mass (Da) Q3 Mass (Da) Dwell(msec) Param Start Stop ID 225.000 42.000 50.00 CE -46.00 -46.00 Pentobarbital 4 CXP -7.00-7.00

@Q1 Mass (Da) Q3 Mass (Da) Dwell(msec) Param Start Stop ID 230.000 187.000 50.00 CE -17.00 -17.00 d5-Pentobarbital CXP -10.00 -10.00

Parameter Table(Period 1 Experiment 1):

CAD: Medium
GS1: 50.00
GS2: 60.00
CUR: 30.00
TEM: 400.00
IS: -3500.00
DP -100.00
EP -10.00

Valco Valve Diverter

FOOD AND DRUG ADMINISTRATION OFFICE OF REGULATORY AFFAIRS Denver Laboratory	Document Number: WI-000140 Compendium method C-005.xx	Revision #: 04 Revised: Nov. 2019
Title: Determination of pentobarbital in ingredients of animal origin and in finished pet foods using liquid chromatography tandem mass spectrometry.		Page 32 of 34

al Time (min)	Position
0.0	В
3.0	Α
5.0	В
	0.0 3.0

Agilent LC Pump Method Properties Pump Model: Agilent 1260 Binary Pump

Minimum Pressure (psi): 0.0 Maximum Pressure (psi): 8702.0

Dead Volume (µI): 40.0

Maximum Flow Ramp (ml/min²): 100.0 Maximum Pressure Ramp (psi/sec): 290.0

Max Flow Ramp Up (ml/min²): 100.0 Max Flow Ramp Dn (ml/min²): 100.0

Step Table:

@Step	Total Time(min)	Flow Rate(µl/min)	A (%)	B (%)
0	0.00	350	95.0	5.0
1	3.50	350	5.0	95.0
2	4.50	350	95.0	5.0
3	8.50	350	95.0	5.0

Left Compressibility: 50.0

Right Compressibility: 115.0 Left Dead Volume (μl): 40.0 Right Dead Volume (μl): 40.0 Left Stroke Volume (μl): -1.0 Right Stroke Volume (μl): -1.0

Left Solvent: A2 Right Solvent: B1

Agilent LC Pump Method Properties Pump Model: Agilent 1260 Binary Pump

Minimum Pressure (psi): 0.0 Maximum Pressure (psi): 8702.0

Dead Volume (µI): 40.0

Maximum Flow Ramp (ml/min²): 100.0 Maximum Pressure Ramp (psi/sec): 290.0 Max Flow Ramp Up (ml/min²): 100.0

FOOD AND DRUG ADMINISTRATION OFFICE OF REGULATORY AFFAIRS Denver Laboratory Title: Determination of pentobarbital in ingredients of animal origin and in finished pet foods using liquid chromatography tandem mass spectrometry. Page 33 of 34

Max Flow Ramp Dn (ml/min²): 100.0

Step Table:

@Step	Total Time	e(min) Flow Rate(µl/min)	A (%)	B (%)
0	0.00	0	50.0	50.0
1	8.50	0	50.0	50.0

Left Compressibility: 50.0

Right Compressibility: 115.0 Left Dead Volume (µI): 40.0 Right Dead Volume (µI): 40.0 Left Stroke Volume (µI): -1.0 Right Stroke Volume (µI): -1.0

Left Solvent: A2 Right Solvent: B2

Agilent Column Oven Properties

Left Temperature (°C): 40.00
Right Temperature (°C): 40.00
Temperature Tolerance +/- (°C): 1.00
Start Acquisition Tolerance +/- (°C): 1.00

Time Table (Not Used)

Column Switching Valve Installed 10Port2Pos

Position for first sample in the batch: Left Use same position for all samples in the batch

CTC PAL Autosampler Method Properties

Loop Volume1 (µI): 20 Loop Volume2 (µI): 20 Injection Volume (µI): 5.000 Barcode Reading: Disabled

Method Description:

Syringe: 100uIDLW

Cycle date: 9/9/2010 3:26:06 PM

Cycle name: Analyst LC-Inj DLW Fast_Rev05

For the most current and official copy, check QMiS.

FOOD AND DRUG ADMINISTRATION OFFICE OF REGULATORY AFFAIRS Denver Laboratory	Document Number: WI-000140 Compendium method C-005.xx	Revision #: 04 Revised: Nov. 2019
Title: Determination of pentobarbital in ingredients of animal origin and in finished pet foods using liquid chromatography tandem mass spectrometry.		Page 34 of 34

Airgap Volume (µI)	3
Front Volume (µI)	5
Rear Volume (µI)	5
Filling Speed (µl/s)	5
Pullup Delay (ms)	3
Inject to I	_C VIv1
Injection Speed (µl/s)	5
Pre Inject Delay (ms)	500
Post Inject Delay (ms)	500
Needle Gap Valve Clean (m	ım) 3
Valve Clean Time Solvent 2	(s) 3
Valve Clean Time Solvent 1	(s) 4
Post Clean Time Solvent 1 ((s) 3